


Non-animal method for predicting skin sensitization (OECD TG442C)

ADRA (**A**mino acid **D**erivative **R**eactivity **A**ssay) is a *in chemico* test method, representing an alternative (non-animal) method for the evaluation of skin sensitization compounds. The kit comprises two amino acid derivatives, NAC (N-(2-(1-naphthyl)acetyl)-L-cysteine) and NAL (α -N-(2-(1-naphthyl)acetyl)-L-lysine), which are composed out of a nucleophilic region and a detection molecule (naphthalene rings).

ADRA is proposed to address the molecular initiating event of skin sensitization AOP (**A**dverse **O**utcome **P**athway) by quantifying the reactivity of the test compounds against the model synthetic amino acid derivatives. The reaction is monitored by HPLC-analysis after 24 hours incubation at 25°C and 281 nm, determining the relative residual concentration of NAC and NAL in the reaction liquid.

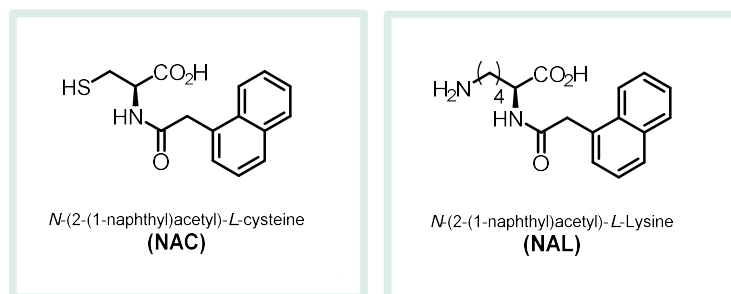

Advantages of ADRA

- NAC and NAL are detected at a relatively long wavelength (281 nm).
 - Preventing co-elution of the test chemical and the nucleophilic reagent.
- ADRA avoids precipitation by using only 1% of reactants of the existing method.
- Variability of result values are very low.
 - In ADRA, the reaction is stopped by the addition of a fixing solution (2.5%TFA) before analyzing
- Multiple test chemicals can be assayed in a short period of time.
 - The test procedure is performed using a single 96-*well* plate and a multichannel pipette.


Kit Reagents



Components	Amount
NAC	2 x 10 ml
NAL	2 x 10 ml
NAC Buffer (pH 8.0) premixed	2 x 300 ml
NAL Buffer (pH 10.2) premixed	2 x 300 ml
0.01 mol/l EDTA Solution	2 x 1 ml

Code No.	Product Name	Package Size
296-80901	ADRA Kit	1 Kit

Additional reagents required

Code No.	Product Name	Package Size
204-02743	Trifluoroacetic Acid 	25 ml
015-08633	Acetonitrile	3 l
217-01031	Ultrapure Water *	1 l
016-00346	Acetone **	500 ml
043-07216	Dimethyl Sulfoxide **	500 ml

* Use water with low metal content.

** Not to be used, if the test chemical dissolves in water or acetonitrile.

Phenylacetaldehyde (CAS RN[®] 122-78-1) is required as a positive control.

Apparatus required

- 1) Electronic balance – precision scale: ± 0.1 mg
- 2) Three Micropipettes – 10 μ l, 100 μ l and 1000 μ l
- 3) 12 channel pipette – 150 μ l
- 4) HPLC system –light-shielding auto-sampler for 96-*well* plates and 0.3 ml/min liquid feeding
- 5) UV detector – photodiode array (PDA) detector or absorbance detector (281 nm)
- 6) HPLC column
- 7) pH meter – precision scale: ± 0.01 pH
- 8) Incubator – 25 °C
- 9) 96-*well* plates
- 10) 500 mL plastic bottles
- 11) Vortex mixer
- 12) Plate seal*
* Use the seal having high sealability and solvent-resistant performance.
- 13) Plate shaker
- 14) Plate centrifuge

Note: to avoid NAC/NAL dimerization upon metal ion contamination, all consumables (except for components for HPLC analysis) must be made of polypropylene or polyethylene.

Listed products are intended for laboratory research use only, and not to be used for drug, food or human use. / Please visit our online catalog to search for other products from FUJIFILM Wako; <https://labchem-wako.fujifilm.com> / This leaflet may contain products that can not be exported to your country due to regulations. / Bulk quote requests for some products are welcomed. Please contact us

FUJIFILM Wako Pure Chemical Corporation

<https://labchem-wako.fujifilm.com>
1-2, Doshomachi 3-Chome Chuo-Ku, Osaka 540-8605, Japan
Tel: +81-6-6203-3741 • Fax: +81-6-6203-1999
ffwk-cservice@fujifilm.com



FUJIFILM Wako Chemicals U.S.A. Corporation

<https://www.wakousa.com/>
1600 Bellwood Road Richmond, VA 23237, USA
Toll-Free (U.S. only): +1-877-714-1920
Tel: +1-804-714-1920 • Fax: +1-804-271-7791
wkuslabchem@fujifilm.com

FUJIFILM Wako Chemicals Europe GmbH

<https://www.wako-chemicals.de>
Fuggerstraße 12, 41468 Neuss, Germany
Tel: +49-2131-311-0 • Fax: +49-2131-311 100
labchem_wkeu@fujifilm.com