

For blood-brain barrier research

Anti Claudin-5, Monoclonal Antibody(R9/M48)

Claudin-5 is a tetra-transmembrane protein with approximately 23kDa, belongs to a member of Claudin family. It functions as a key component regulating tight junction of blood-brain barrier¹⁾. These products are monoclonal antibodies which specifically recognize Claudin-5, and have neutralizing activity which reduces barrier function in blood-brain barrier cell model¹⁾. They have no cross-reaction to Claudin-1-4,6,7¹⁾.

1) Hashimoto, Y., et al.: *J. Pharmacol. Exp. Ther.*, 363(2), 275(2017).

Antibody information

	Anti Claudin-5, Monoclonal Antibody (R9)	Anti Claudin-5, Monoclonal Antibody (M48)
Code No.	014-28101	011-28091
Clone No.	R9	M48
Antigen	Vector expressing human claudin-5 full length accession#AAH19290	Vector expressing human claudin-5 full length accession#AAH19290
Buffer	PBS	PBS
Subclass	Rat IgG2b · κ	Mouse IgG3 · κ
Cross-reactivity	Human, monkey	Human, monkey
Specificity	Claudin-5 No cross-reaction to Claudin-1-4,6,7 ¹⁾	Claudin-5 No cross-reaction to Claudin-1-4,6,7 ¹⁾
Concentration	Indicated to the label First lot : 1.1mg/mL	Indicated to the label First lot:1.2mg/mL
Application	Flow cytometry 0.5-5µg/mL Neutralizing assay(in vitro) 90-150µg/mL	Flow cytometry 0.5-5µg/mL Neutralizing assay(in vitro) 90-150µg/mL


Product Name	Wako Cat. No	Pkg. Size	Grade	Storage Condition
Anti Claudin-5, Rat Monoclonal Antibody (R9)	014-28101	500µL	For Immunochemistry	Keep at -80°C
Anti Claudin-5, Monoclonal Antibody (M48)	011-28091	500µL	For Immunochemistry	Keep at -80°C

This product is sold under the license of Osaka University-Ehime University-FUJIFILM Wako Pure Chemical Corporation joint application patent. PCT/JP2018/016975 「Anti-cldn-5 antibody, and drug containing said antibody」

Application data (Neutralizing assay *in vitro* BBB cell model)


The barrier activity was evaluated by adding these antibodies (R9 and M48) were added to the blood-brain barrier cynomolgus cell model (BBB kit, PharmaCo-cell).

(1) Trans-epithelial/endothelial electrical resistance (TEER)

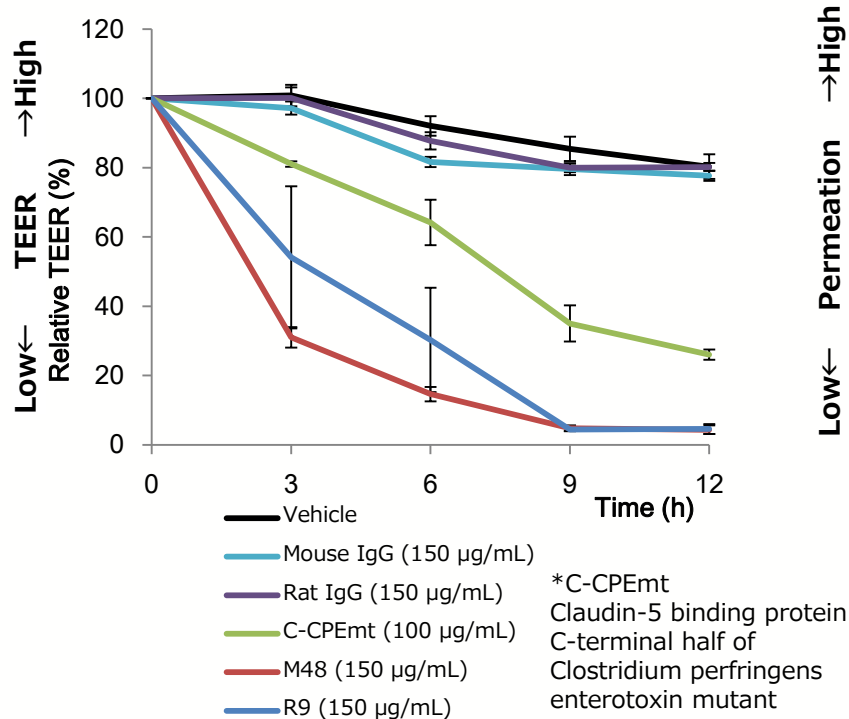

These antibodies (R9 and M48) reduced TEER. The effects were higher than C-CPEmt.

(2) Substance permanent

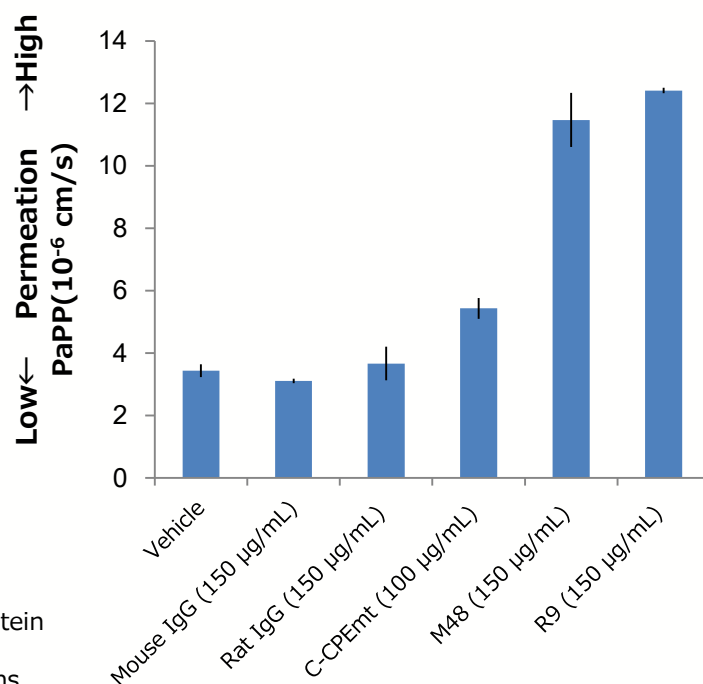

These antibodies (R9 and M48) increased permanent of fluorescein isothiocyanate-labeled dextran (4kDa). The effects were higher than C-CPEmt.

Listed products are intended for laboratory research use only, and not to be used for drug, food or human use. / Please visit FUJIFILM Wako Laboratory Chemicals site: <https://labchem-wako.fujifilm.com/> / This leaflet may contain products that cannot be exported to your country due to regulations. / Bulk quote requests for some products are welcomed. Please contact us.

FUJIFILM Wako Laboratory Chemicals site
<https://labchem-wako.fujifilm.com>



FUJIFILM Wako Pure Chemical Corporation
 1-2, Doshomachi 3-Chome, Chuo-ku, Osaka 540-8605, Japan
 Tel: +81 6 6203 3741 Fax: +81 6 6203 1999
 ffwk-cserservice@fujifilm.com

FUJIFILM Wako Chemicals U.S.A. Corporation
 1600 Bellwood Road, Richmond, VA 23237, U.S.A.
 Toll-Free (U.S. only): +1 877 714 1920
 Tel: +1 804 271 7677 Fax: +1 804 271 7791
 wkuslabchem@fujifilm.com

FUJIFILM Wako Chemicals Europe GmbH
 Fuggerstr 12, 41468 Neuss, Germany
 Tel: +49 2131 311 0 Fax: +49 2131 311 100
 labchem_wkeu@fujifilm.com